



A SIGN OF QUALITY


## HOW TO DOWNLOAD DATA LOGGER ON NLMB1006

1. Insert USB stick into front panel.



2. Press and hold down arrow and light bulb. USB with display on the screen, Keep holding until USB is no longer on the screen and the temperature is displayed. Remove USB stick

3. Open File on computer

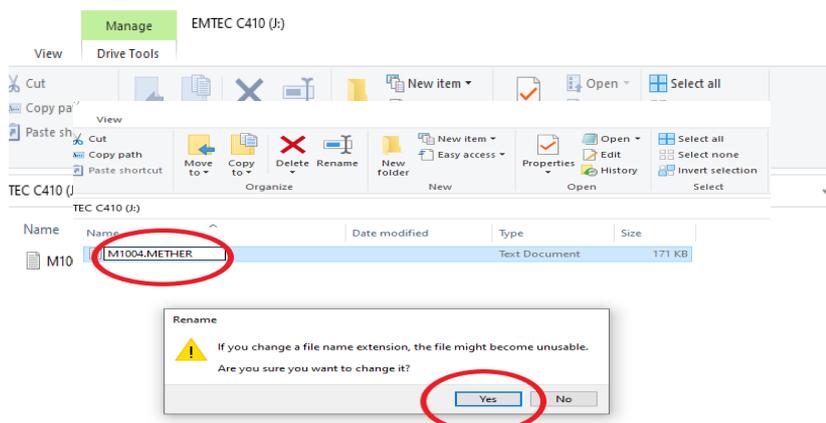

4. Change file name to **.METHER** – then press **Yes**




A SIGN OF QUALITY

5. Open Software



6. 1- Change to English
- 2- Choose Single Temperature
- 3- Choose whether you want to save as PDF or Excel
- 4- Press Import
- 5- Find **METHER** in D Drive
- 6- Click on file and open
- 7- Data will import then click Okay
- 8- Save AS (Where you want to save)

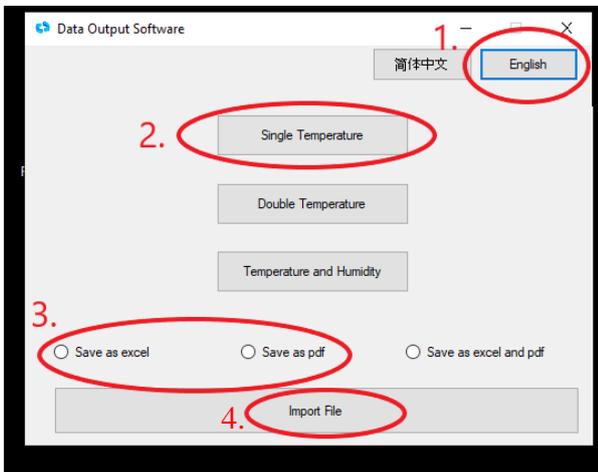



A SIGN OF QUALITY

5. Find where you saved file and it should open up like this.

	A	B	C	D	E
778	2021/12/07	20:36	05.4		
779	2021/12/07	20:46	05.6		
780	2021/12/07	20:56	05.4		
781	2021/12/07	21:06	05.6		
782	2021/12/07	21:16	05.3		
783	2021/12/07	21:26	05.6		
784	2021/12/07	21:36	05.4		
785	2021/12/07	21:46	05.8		
786	2021/12/07	21:56	05.3		
787	2021/12/07	22:06	05.8		
788	2021/12/07	22:16	05.4		
789	2021/12/07	22:26	05.6		
790	2021/12/07	22:36	05.4		
791	2021/12/07	22:46	05.6		
792	2021/12/07	22:56	05.4		
793	2021/12/07	23:06	05.8		
794	2021/12/07	23:16	05.6		
795	2021/12/07	23:26	05.6		
796	2021/12/07	23:36	05.6		
797	2021/12/07	23:46	05.5		
798	2021/12/07	23:56	05.8		
799	2021/12/08	00:06	05.5		
800	2021/12/08	00:16	05.6		

Distributor  
**EVOCARE AUSTRALIA PTY LIMITED**

A.B.N. 98 078 566 604

Trading as **EVOCARE** and **L&M EQUIPMENT**

P.O. Box 145, Everton Park Qld. 4053

Ph: 07 3355 8000

Website: <http://www.evocare.com.au>

Email: [sales@evocare.com.au](mailto:sales@evocare.com.au)

[workshop@evocare.com.au](mailto:workshop@evocare.com.au)

[warehouse@evocare.com.au](mailto:warehouse@evocare.com.au)

[accounts@evocare.com.au](mailto:accounts@evocare.com.au)

